


Quick Guide

Monkey NGAL ELISA Kit (KIT 045)

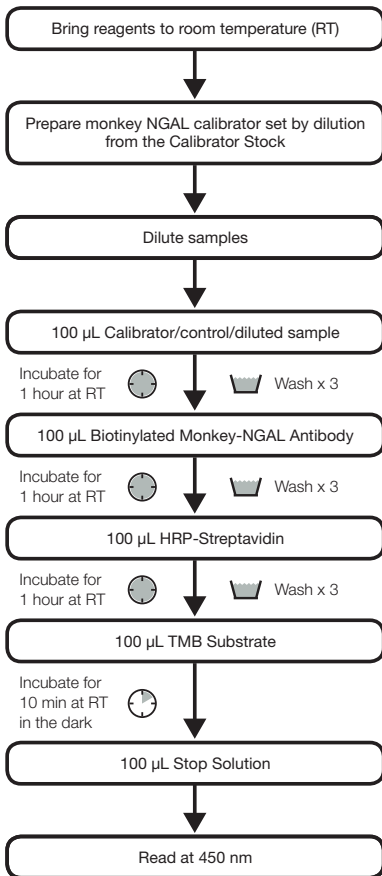


BIOPORTO[®]
Diagnostics


CALBRATOR DILUTION GUIDE*

Monkey NGAL Calibrator Stock

5 mL 1x sample diluent

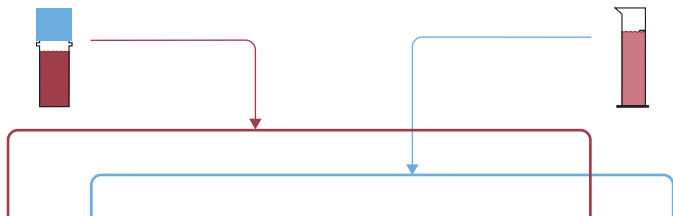

Dilution Factor	1	2	4	10	20	40	100	-
Monkey NGAL	200 µL	200 µL	100 µL	40 µL	20 µL	10 µL	10 µL	-
Sample Diluent	-	200 µL	300 µL	360 µL	380 µL	390 µL	990 µL	400 µL

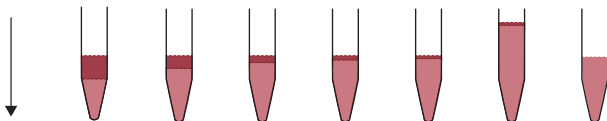

Calibrator #	8	7	6	5	4	3	2	1
NGAL Concentration	1000 pg/mL	500 pg/mL	250 pg/mL	100 pg/mL	50 pg/mL	25 pg/mL	10 pg/mL	0 pg/mL

Transfer to wells	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2	100 µL/ well x 2
-------------------	---------------------	---------------------	---------------------	---------------------	---------------------	---------------------	---------------------	---------------------

*Calibrators are to be freshly diluted immediately prior to use.